


11. TROUBLE SHOOTING

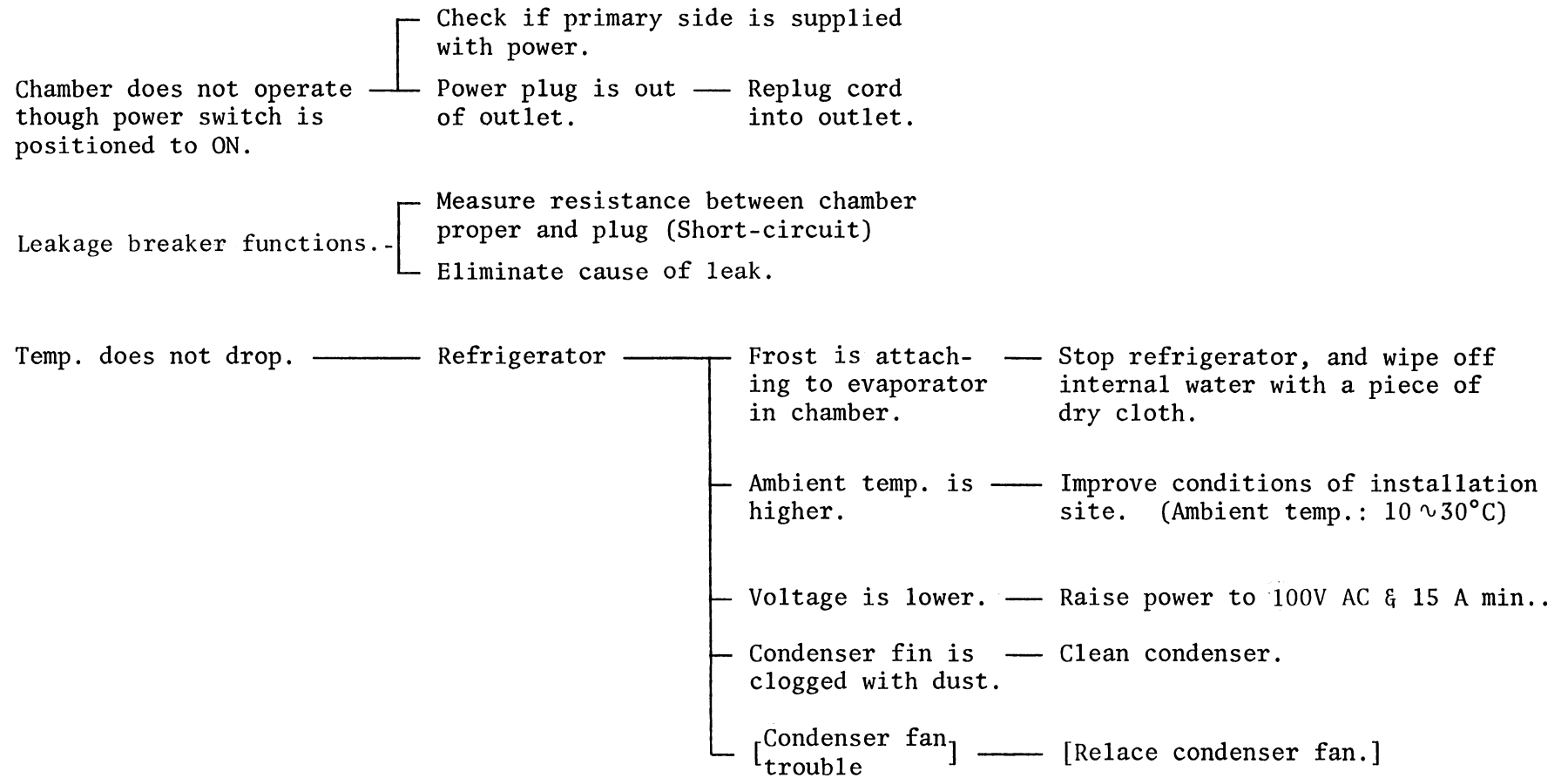

Temp. control is not done.

In spite of turning temp. controller knob, heating heater pilot lamp keeps light or keeps dark.

Thermal fuse sensor is broken.

Replace thermal fuse sensor.

Power voltage is lower.

Raise power to 100V AC & 15 A min..

Ambient temp. is higher.

Improve conditions at installation site.

Automatic temp. over-heat protection circuit functions.

Reset.
(Set 5~10°C higher than setting level of controller.)

Air circulator is turning.

Check setting temp..

◦Digital display section is blinking.

Temp. indication blinks at 999°C.

Repalce temp. sensor.

Remote input voltage is not adequate.

Temp. indication blinks at -70°C and under.

In spite of remote operation, remote input/output terminal does not receive signal.

In case of local operation, position remote SELECT switch to LOCAL.